



Cod UNG in RT-LAMP

Carry-over
contamination control

www.arcticzymes.com

Removal of carry-over contamination in RT-LAMP

ArcticZymes Cod UNG is ideal for contamination control in RT-LAMP. One unit of Cod UNG per 30 µl reaction is sufficient for removing even high concentrations of carry-over contamination.

- ✓ Ensure that dNTP mixes containing dUTP were used in previous RT-LAMP experiments.
- ✓ Check that the RT-LAMP reaction is compatible with dUTP by running side-by-side reactions containing different ratios of dUTP to dUTP (100% dUTP, 90% dUTP, 80% dUTP and 0% dUTP).
- ✓ Add 1 U Cod UNG per 30 µl reaction in the RT-LAMP reaction mix, see suggested RT-LAMP reaction below. Prepare reaction mix on ice.
- ✓ Analyze your RNA at 65°C, no preincubation is necessary.

Additional information

ArcticZymes Cod UNG is compatible with post-amplification analysis as it is completely and irreversibly inactivated at 55°C.

Suggested RT-LAMP reaction

Component	1 rxn (µl)	Final Concentration
IsoPol™ BST+ 10X Reaction Buffer Pack	3	1X
dNTP mix (10 mM, dTTP replaced with dUTP)	4.2	1.4 mM each
MgSO ₄ (100 mM)	1.2	4 mM (8 mM total) *
EvaGreen® Fluorescent DNA Stain, 100 µM	0.39	1.3 µM
IsoPol™ BST+, 8 U/µl	1.2	0.32 U/µl
FIP + BIP 15 µM/each F3 + B3 3.75 µM/each LF + LB 7.5 µM/each	1.6	0.8 - 1.6 µM 0.2 µM 0.4 µM
AMV-RT (21 U/µl)	0.09	0.063 U/µl
ArcticZymes Cod UNG (5 U/µl)	0.2	1U (0.033 U/µl)
PCR grade water	13.12	N/A
RNA template	5	N/A
Final volume	30	

*4mM contribution from the IsoPol™ BST+ 10X Reaction Buffer

Disclaimer: These products are intended for further manufacturing use or research use only. Certain applications of ArcticZymes Technologies ASA products may require licenses from others. It is the expressed duty of any receiver of ArcticZymes Technologies ASA products to acquire such licenses, if necessary. To the extent allowed by law, ArcticZymes Technologies ASA will not be liable for damages, whether direct, indirect, incidental, or consequential in connection with or arising from this document, including the use of it. ArcticZymes Technologies ASA products may be covered by pending or issued patents, designs or design applications and/or trademarks or trademark applications or any other registered or unregistered Intellectual Property Right. Version 1.0 • October 2021

ArcticZymes Technologies ASA

Sykehusveien 23
N-9294 Tromsø, Norway

T (47) 7764 8900
E contact@arcticzymes.com
I www.arcticzymes.com

ArcticZymes Technologies ASA

987 Old Eagle School Road, Suite 709
Wayne, PA 19087, USA

T (484) 534 3567
F (484) 368 3558
E contact-us@arcticzymes.com

ArcticZymes AS

Handelswei 1A 8501 XJ
Joure, the Netherlands

T (47) 7764 8900
E contact@arcticzymes.com
I www.arcticzymes.com

